


BRO/BUF 08

A laboratory pilot plant designed for process development, membranes scale up testing, quality assurance and small-scale production.

A simple self-contained unit of minimum hold-up volume (15 Liters) with electric motor, CAT 1051 high pressure pump, membrane module, heat exchanger, pressure gauges and flow meter. All mounted on a welded stainless-steel framework.

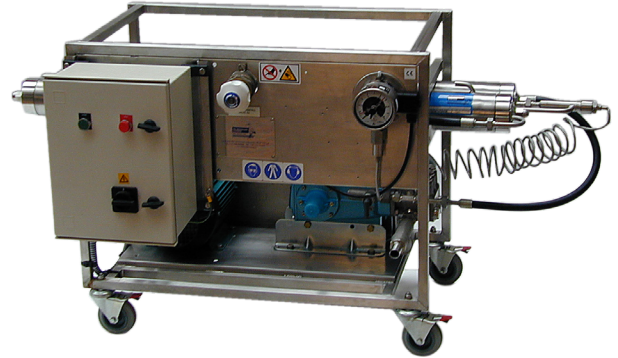

Application Range

Microfiltration (MF)	Ultrafiltration (UF)	Nanofiltration (NF)	Reverse Osmosis (RO)
up to 30 L/min	up to 30 L/min	18 – 22 L/min	18 – 22 L/min

Electrical Power Supply	Connections
400 Volts	Feed inlet to pump: 3/4" OD hose
3-phase (5 pins)	HE in/out cooling water: 1/2" OD hose
50 Hz. 32 Amp	Permeate & Shroud outlet: 1/2" OD hose
Motor rating: 5.5 KW	Concentrate outlet UF: 3/4" hose
	Concentrate outlet RO: 1/2" hose

Size & Weight

Dimension			Unit & Package	Unit Only
Length	Height	Width		
1,9 m	1,05 m	0,8 m	350 kg	244 kg

Framework

Welded stainless-steel frame fabricated from a high grade SS (AISI 304) for corrosion resistance, rigidity and lightness.

Module

4 ft (1.2 m) B1 Module

Comprised of 18 perforated stainless-steel tubes in the form of a shell & tube, each tube fitted with a membrane element. The module is designed with a series flow end caps; connecting all 18 tubes in series.

- Membrane area: 0.88 m²
- Module weight: 14.4 kg
- Hold-up volume: Tubeside 2.8 L, Shroud-side 6.7 L
- Membrane tube ID: 12.7 mm

Heat Exchanger (HE)

Shell & Tube Type Heat Exchanger 2 Ft (0.6 M)

Process fluid is piped from the outlet of the pump through all 18 tubes in series within the heat exchanger while cooling water passes at low pressure (1-3 bar) through the shroud (shell) side. Cooling media flow could be up to 20 L/min.

Module tube side mechanical operating limit

- Max operating pressure: up to 64 bar
- Max pressure drop: 10 bar
- Max operating temperature: up to 80°C

2020 11 11 DS

Corporate Office

PCI Membranes
Unit 11, Fulcrum 2, Solent Way, Whiteley, Fareham,
Hampshire, PO15 7FN, United Kingdom

Phone

Phone: + 44 1489 563 470
Email: pcimembranes@filtrationgroup.com


Online

Website: www.pcimembranes.com
www.filtrationgroup.com

BRO/BUF 08

A laboratory pilot plant designed for process development, membranes scale up testing, quality assurance and small-scale production.

A simple self-contained unit of minimum hold-up volume (15 Liters) with electric motor, CAT 1051 high pressure pump, membrane module, heat exchanger, pressure gauges and flow meter. All mounted on a welded stainless-steel framework.

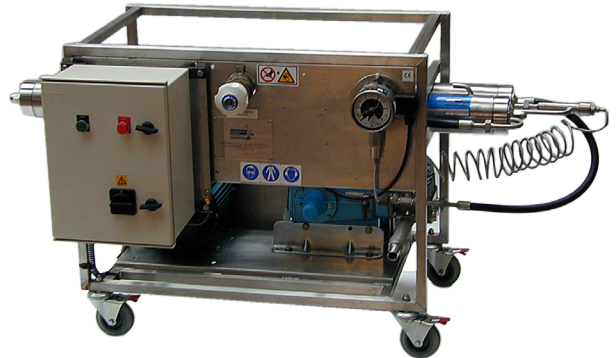

Pump

- A CAT 1051 HP (piston type) is fitted.
- Capable of a flow rate between 6 – 38 L/min.
- Flow can be adjusted via a VSD on the panel.

Motor & Drive Guard

- 5.5 KW motor with 3 phase, 1 earth and 1 neutral (5 pins).
- TEFC foot mounted motor with simple adjustment for drive belt tensioning and alignment. A cover for the toothed belt drive between pump and motor is also fitted.

Pressure Relief Valve

- The plant can either be fitted with a pressure relief valve set at 70 bar for NF/RO or 20 bar for MF/UF operation.

Pulsation Damper

A pulsation damper or accumulator is fitted to the pump outlet. Charge with Nitrogen as follows:

- 40 bar for RO.
- 6 bar or below for MF/UF operation (depending on the operating pressure).

Strainer

- A 40-mesh (400 micron) “Y” type strainer is fitted to the pipe work to protect the pump and modules.

Flowmeter

- Variable Area PN40 flowmeter is fitted on the feed line.

Pressure & Temperature Sensor. Pressure Control Valve

- Both module inlet and outlet pressure are constantly monitor and display on independent pressure gauges.
- A hand operated needle valve for NF/RO operation or a diaphragm valve for MF/UF operation is used to create the back pressure.

Safety

- The unit is designed to the principles of Supply of Machinery (Safety) Regulations 1992 and are safe if operated in accordance with the procedures in the operating manual.

Corporate Office

PCI Membranes
Unit 11, Fulcrum 2, Solent Way, Whiteley, Fareham,
Hampshire, PO15 7FN, United Kingdom

Phone

Phone: + 44 1489 563 470
Email: pcimembranes@filtrationgroup.com

Online

Website: www.pcimembranes.com
www.filtrationgroup.com